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🇬🇧 Cleaning & Maintenance

1. Clean b.safe Waste Caps with a damp cloth and, if necessary, with detergents and warm water. **Attention:** Water must not get into the b.safe Exhaust Filter. We recommend to remove it before cleaning the Cap.
2. Possible labelling or adhesive residues on b.safe Waste Caps can be removed with isopropyl.
3. Once the b.safe Service Life Indicator is activated, the white inspection window turns into red with increasing time of use. If the white writing "CHANGE" is completely visible on the red background, the b.safe Exhaust Filter is dead and has to be replaced.

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b.safe Abluftfilter / b.safe Exhaust Filter

| Größe Size | Standzeit Service life | Verpackungseinheit Packing unit | Artikelnummer Cat. No. |
|---------------|---------------------------|------------------------------------|---------------------------|
| Small | 3 Monate / months | 2 Stück / pieces | M 506-01 |
| Medium | 6 Monate / months | 2 Stück / pieces | M 506-02 |
| Large | 6 Monate / months | 2 Stück / pieces | M 506-03 |

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Push the capillaries through a suitable b.safe Fitting and afterwards through the b.safe Waste Cap.

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Push the capillaries through the b.safe Fittings until their ends protrude on the bottom side of the b.safe Waste Cap for about 10 mm. Tighten the b.safe Fittings.

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Screw the b.safe Hose Connector into the suitable thread of the b.safe Waste Cap. Push the tubing on the hose connector until it sticks.

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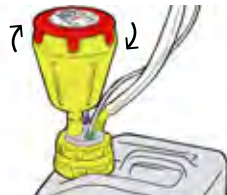
Tighten the b.safe Waste Cap onto a suitable canister until the insert does not move freely any longer.

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Screw a b.safe Exhaust Filter into the free thread connection of the b.safe Waste Cap.

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Push the button  to activate the Service Life Indicator. Optional: Note the exchange date on the label.

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Mark the b.safe Waste Cap with a foil pen or an adhesive label if necessary.

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Important: Use b.safe Blind Fittings for gas tight closing of unused openings on a b.safe Waste Cap!

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Bohlender GmbH · Waltersberg 8 · 97947 Grünsfeld